

HIGH VOLUME SLIDERITE®


SECTION A

GENERAL USE

The Sliderite® is a mobile pallet developed to assure that the Cryoport High Volume shipper's remains in the upright position.

Maintaining this orientation assures that the shippers validated temperature hold time can be achieved.

It also allows for easy mobility of the High Volume Shipper.

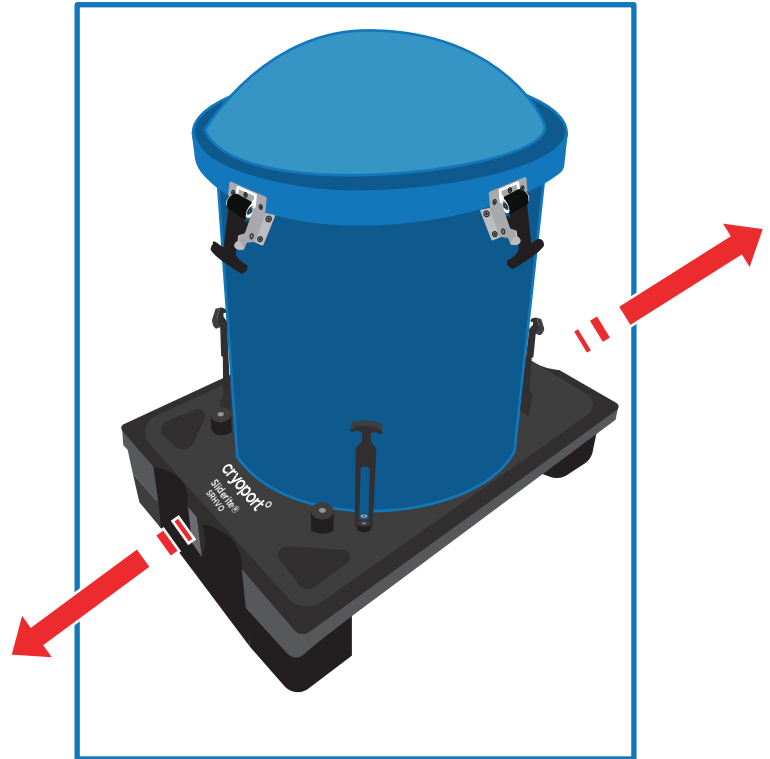


SECTION B

MODES

The Sliderite® pallet has two primary modes that are controlled by the pedals on either end of the pallet.

Pushing down on the **RED** pedal puts the pallet into the stationary mode. In this mode the pallet will not roll.

Pushing on the **GREEN** pedal will release the pallet into the motion mode and allows the pallet to be rolled around.

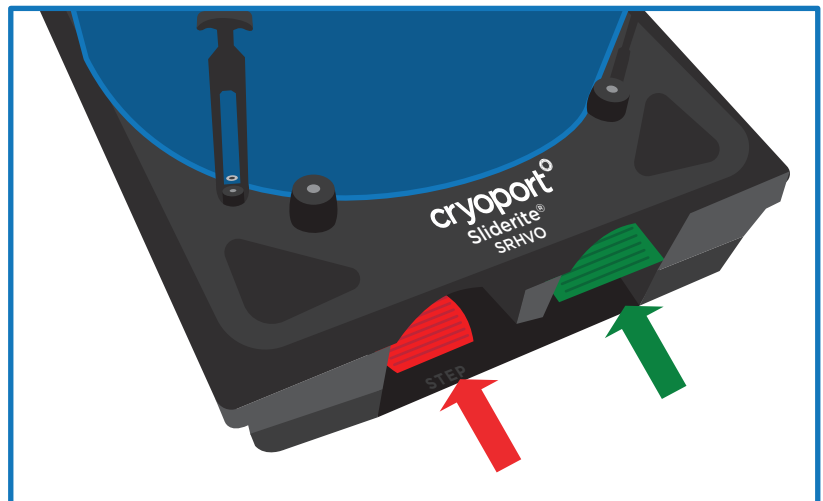

IF THE DRY SHIPPER DOES NOT EMIT VAPOR WHEN PLUG IS REMOVED

PLEASE CONTACT CUSTOMER SERVICE IMMEDIATELY, BY CALLING (949) 470-2305 OR EMAIL CS@CRYOPORT.COM

24-hour Help Line
(949) 470-2305

Cryoport, Inc. • 17305 Daimler St. • Irvine, CA 92614
Phone: 1-949-470-2300 | Fax: 1-949-470-2306
www.cryoport.com • e-mail contact: cs@cryoport.com