

### How to Access Digital Data Logger—CXHVO

**Standard Shipment: CXHVO**  
**Digital Hobo Data Logger Access**

**Data Station instructions and software can be found at:**  
<http://www.cryoport.com/data-station-directions>.

**Note:** The data logger will not record internal temperature while disconnected from the thermocouple.

#### Step 1



Open Shipper.

#### Step 2a



Unlatch data logger box using thumbs to press against the black hinge.

#### Step 2b




Open the lid and expose the data logger in its metalized pouch.

#### Step 4



To view temperature data remove the datalogger from the metalized bag. If you wish to download data go to step 5a. If not go to step 5b

#### Step 5a

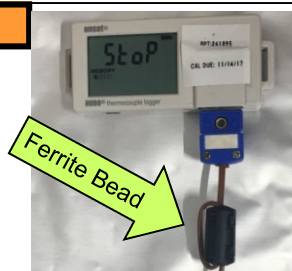


Follow the Data Station instruction to download temperature data, if desired. The data logger may be detached from the thermocouple to move to a remote computer to download data. Once completed go to step 5b

#### Step 5b



Place data logger back into the pouch and make sure to push the cable with the ferrite bead completely down into the bag. Push the cable into the far corner of the bag and reseal by pushing all the way across the zip lock seal.

#### Step 6

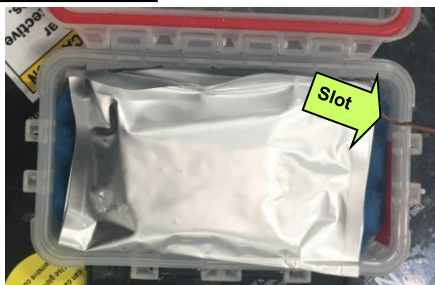

Fold the pouch in half and put the unit back into the protective box. Ensure that the cable is inserted into the slot to prevent cable pinching

#### Step 7



Close and latch the lid.

#### Step 8

Follow the unloading instructions to complete the receiving process and return the shipper to Cryoport.

**cryoport**  
 Standard Shipment: Unloading

**Sample Receiver**  
 How to Unload Samples

**Step 1:** Tear backflaps on Tyvek bags. Remove bag completely.

**Step 2:** Remove sample container.

**Step 3:** Reinsert shipper plug. Push plug fully into place.

**Step 4:** If any slot was used in this area for your shipment please place the enclosed "EMPTY" label over it.

**Step 5:** Remove "STOP" insert from inside shipping envelope on flap "B". Then empty ONLY the top insert!

**24-HOUR HELP LINE**  
 (949) 470-2306

**STOP**

**Instructions Contained on Back of Sheet**

Cryoport, Inc. • 17305 Daimler St. • Irvine, CA 92614  
 Phone: 1-949-470-2300 Fax: 1-949-470-2306  
 www.cryoport.com • E-Mail: cs@cryoport.com

**24-HOUR HELP LINE**  
**(949) 470-2305**

**Cryoport, Inc. • 17305 Daimler St. • Irvine, CA 92614**  
 Phone: 1-949-470-2300 Fax: 1-949-470-2306  
 www.cryoport.com • E-mail Contact: cs@cryoport.com