

### How to Access Hobo Data Logger CXHVO


#### Standard Shipment: CXHVO Standard Hobo Data Logger Access

#### Step 1



Open Shipper.

#### Step 2a

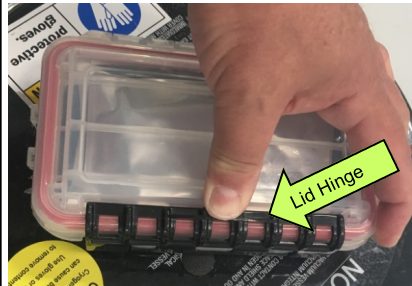

Unlatch Data Logger box using thumbs to press against the black hinge.

#### Step 2b



Open the lid and expose the Data Logger in its metalized pouch.

#### Step 4



Remove the Data Logger from the metalized bag. If you are not downloading data, Go to Step 5a. If your Data Station Software is installed on a nearby computer, Go to Step 5a. Otherwise Go to Step 5b.

#### Step 5a




You can now follow the Data Station instruction to download temperature data, if needed. If Data Station Software is not installed on a nearby computer, disconnect Thermocouple from Data Logger. You can now follow the Data Station instruction to download temperature data.

#### Step 5b



Place Data Logger back into the pouch and make sure to push the cable with the ferrite bead completely down into the bag and push the cable into the far corner of the bag and reseal by pushing all the way across the zip lock seal.

#### Step 6

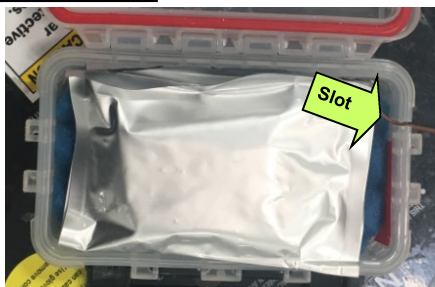

Fold the pouch in half and put the unit back into the protective box. Assure that the cable is inserted into the slot to prevent cable pinching

#### Step 7



Close and latch the lid.

#### Step 8

Follow the Unloading Instructions to complete the receiving process and return the shipper to Cryoport.

**24-HOUR HELP LINE**  
(949) 470-2306

Cryoport, Inc. • 17305 Daimler St. • Irvine, CA 92614  
Phone: 1-949-470-2300 Fax: 1-949-470-2306  
www.cryoport.com • E-mail Contact: cs@cryoport.com