

1. Email Data

Internal Temperature with Standard Deviation may be obtained directly from the email that was provided during Pickup/Delivery in the Details Section or by selection any [Live View Link](#).

Pick Up/Delivery Details Leg X (Company to Company) on order XXXXX has been picked up or delivered and is now en route.

Time of Pickup: XX/XX/XX; X-XX AM/PM -XXXX

Temperature statistics while waiting pickup on leg X, from 11/29/2016 08:48 AM(UTC -0600) to 11/29/2016 03:06 PM(UTC -0600).

Package	Data Logger	Metric	Min	Max	Average	Standard Deviation	Last Reading	Last Reading At	Live View
1	99000512004539	Internal Temperature	-194.75 °C	-194.05°C	-194.38°C	0.18°C	-194.45°C	11/29/2016 1:05 PM (UTC -0600)	Live View


The "Live View" feature may be viewed by clicking on the link provided.
Link will only be viewable if "SRV-LIV" was purchased!

2. Live View

ONLY AVAILABLE IF THE LIVE VIEW SRV-LIV WAS PURCHASED

Important Live View Indicators

Internal Temperature Orientation Battery Level Location

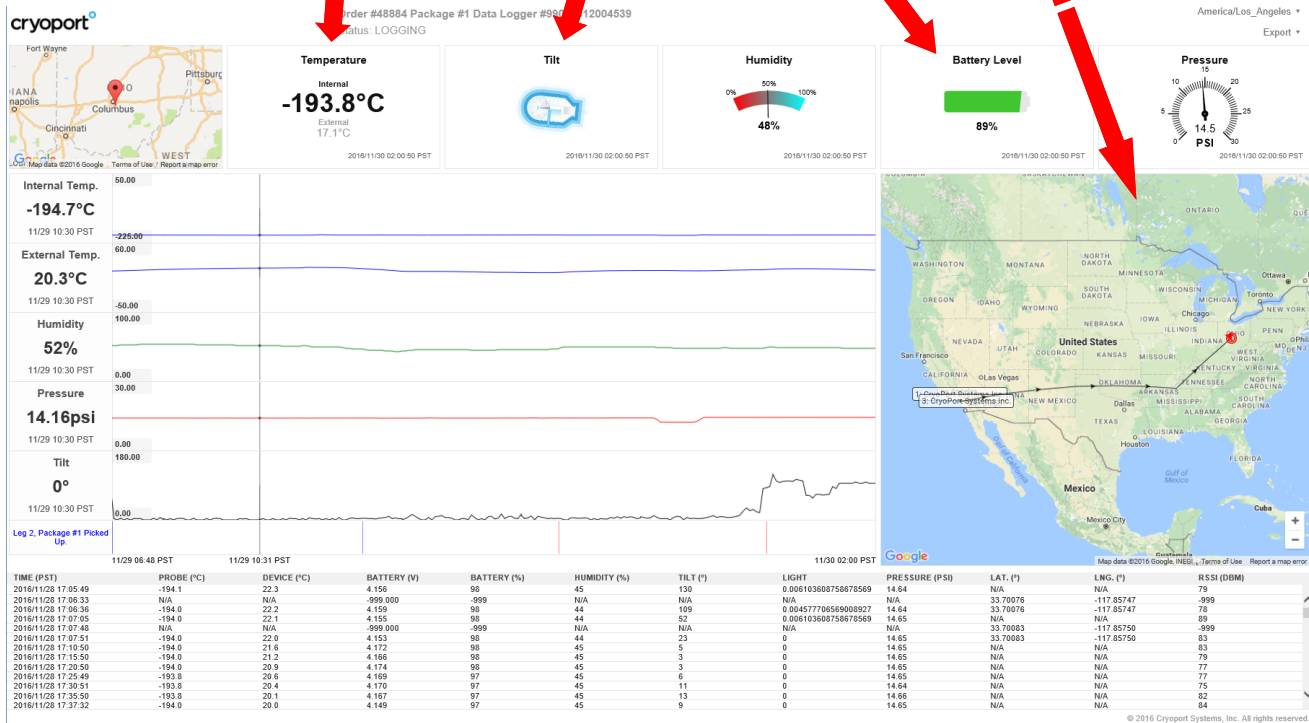

**24-HOUR HELP LINE
(949) 470-2305**

**Instructions Continued
on Back of Sheet.**

3. Portal

Link will only be viewable if "SRV-LIV" was purchased!

The "Live View" feature may be viewed from the portal by searching for the Order Number. Follow these instructions!

B. Search for Order Number

A. Select Order History

C. Select Temperature Monitoring

Order # 49656 for Kathryn Harper Jezwinski (KATH051)
 Phone: 501-454-2758 Fax: Address: 306 Bunnell Street Anchorage AK, 99508 US
 Sales Representative: IVF/Reproductive Medicine Submitted At: 12/13/2016 10:54 AM (UTC-0800)
 Cryoport PO#: Jezwinski_Donor #13964
 Quote #: 13592
 Status: Pending
 Ship From: SEATTLE REPRODUCTIVE MEDICINE, 1505 WESTLAKE AVE N STE 400, SEATTLE WA, 98109-6211 US
 Contact: Jobb, Melanie - 206-301-5015 - melanie.jobb@integramed.com

D. Select Line Item

E. Select Live View

F. Data May be Exported to Excel by selecting Export Data

Package #	Content	Data Logger S...	Dewar Serial ...	Battery Voltage	Sampling Inter...	Delayed Start	Location Interval	Transmit Rate	Next Transmit	Status	Holding Status	Manual Upload
1	Cryopreserve...	99000512004...	24675	4.088	5	No Delay	180	180	11 minutes	Live Logging	PENDING	No

G. Live View will be displayed, see section 2 on Page 1

**24-HOUR HELP LINE
(949) 470-2305**

**Cryoport, Inc. • 17305 Daimler St. • Irvine, CA 92614
 Phone: 1-949-470-2300 Fax: 1-949-470-2306
 www.cryoport.com • E-mail: cs@cryoport.com**