

CRYOPORT ELITE™ 28L ULTRA COLD ADVANCED THERAPY SHIPPER

How to Unload Samples for Shipment

PLEASE FOLLOW LOCAL RULES OF SAFE HANDLING AND DISPOSING OF DRY ICE

Step 1

Wear the appropriate personal protective equipment (PPE) for handling extreme cold temperatures for the unloading of the payload and dry ice.



Step 2



Open the document enclosure lid by pressing down on the lid while simultaneously pressing in each latch on the side one at a time, and then lifting the lid up.



Once document enclosure lid is open, remove blue lid wedge and close the lid by pressing down one side at a time.

Step 3



Prior to removal of serialized zip tie, record serialized number onto all pertinent documentation. Remove serialized zip tie with wire/cable cutters.

Step 4




The lid must be opened one side at a time. Using one hand, press down on the lid while simultaneously pressing the latch release button. Pressing down on the lid will reduce resistance on the button and more easily allow the latch to release.

Step 5



Once lid is open, utilize lid wedge to keep lid propped open and prevent personal injury. The lid wedge should be placed by sliding inward from either side position of the hinge (located between the lid and the base). Lid wedge is located in the document enclosure.

Step 6

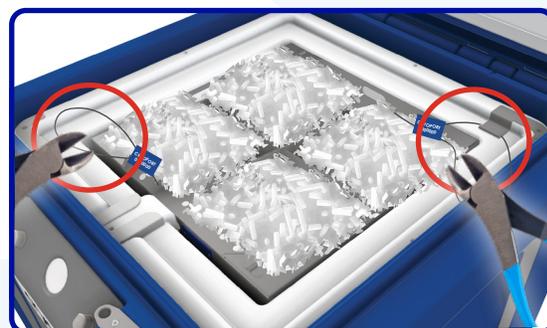

Prior to removal of serialized zip ties, record serialized number onto all pertinent documentation. Remove serialized zip tie with wire/cable cutters.

